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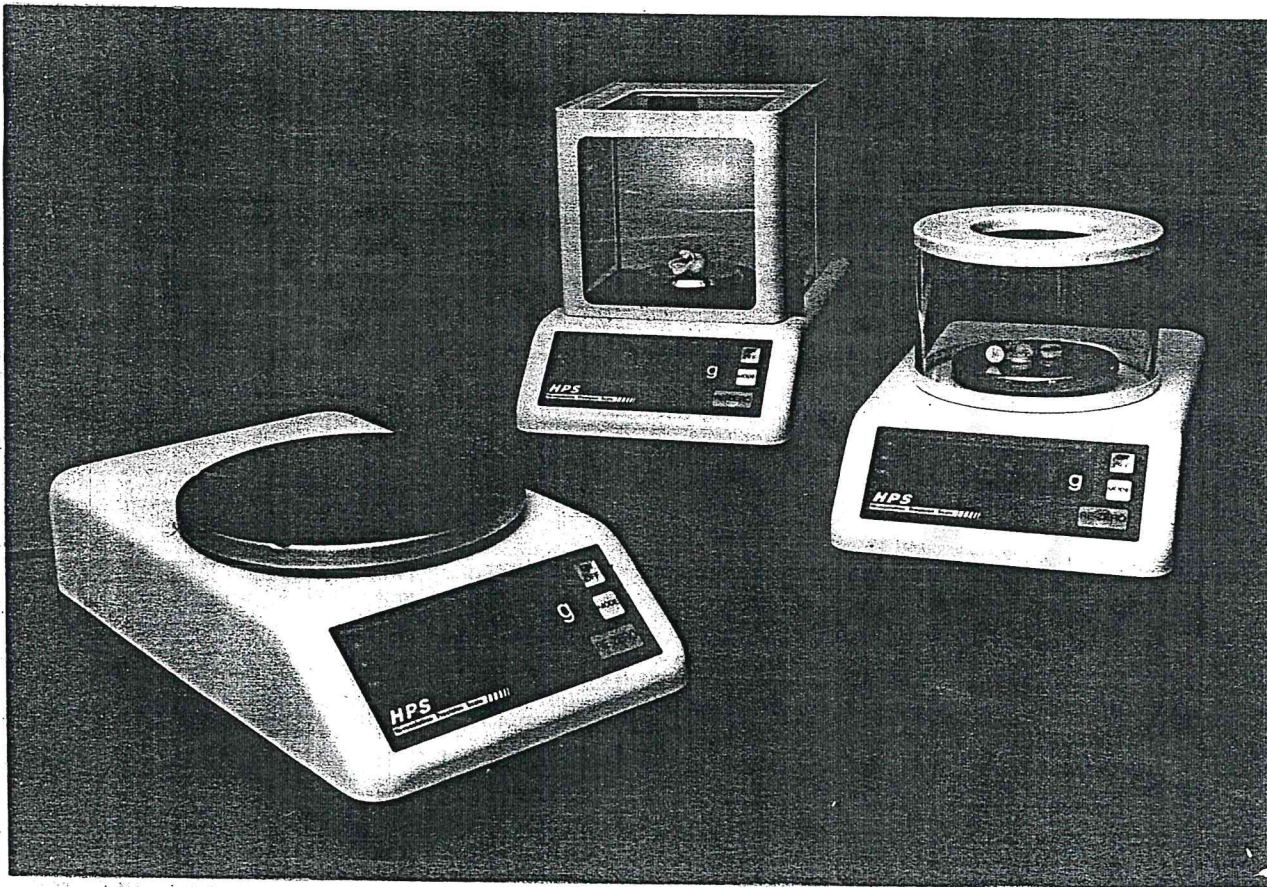
WARNING

N.B.: This "WARNING" is an integral part of the instrument.

This instrument is to be used STRICTLY for scientific purposes and/or for internal factory control ONLY.

Not for the uses referred to in Art. 1 (2) (a) - L. 90/384 EEC.

THIS INSTRUMENT IS NOT TO BE USED IN AREAS WHERE A DANGER OF EXPLOSION EXISTS.

It is important to note the following points for correct operation of the instrument:

- 1 Read the technical instructions given in the operating manual carefully. This manual contains information about installation and use. If these instructions are not carried out carefully, the accuracy of results may be impaired. If in doubt, please consult the supplier's engineers.
- 2 As for any other electrical equipment, installation should be carried out by a competent person. In particular:
 - ensure that the instrument is correctly earthed;
 - do not install the instrument in areas of high fire risk, for example, in the presence of inflammable gases and vapours;
 - switch off and disconnect the power supply before removing any cover;
 - do not touch the instrument with wet hands.
- 3 Install the instrument on a vibration free base and away from draughts or sources of heat. Check the instrument with test weighings after installation or after repositioning.
- 4 Avoid connection to a power supply showing high voltage variations due to other loads. The instrument should not be connected to the same electrical circuit as other high power consuming equipment as large fluctuations may cause some inaccuracy.
- 5 The instrument should be checked every 6 months for correct functioning calibration and accuracy of the internal calibration mass by a suitably qualified person.
- 6 A routine calibration check should be carried out each day before use, using a standard mass.
- 7 Check the zero function before each weighing operation.
- 8 If in doubt about any possible malfunction please call a qualified person. Avoid interference by non-qualified people.
- 9 If the instrument is used for weighing food or drink, ensure that other substances are not weighed on the unit which might cause contamination.
- 10 If the instrument is used for weighing chemicals and other similar substances ensure that it is cleaned and any spillage removed to avoid potentially dangerous reactions.
- 11 Always clean the instrument after use to avoid subsequent inaccuracies and/or damage to the equipment.
- 12 s.r.l. does not accept any liability for damages or other consequences due to tampering of the instrument or its accessories.

ADDITIONAL PROCEDURES FOR WEIGHING SUBSTANCE WITH GREAT ACCURACY

Instruments are manufactured to be reliable in use with high accuracy. However, reduced accuracy is always possible due to ambient conditions, incorrect installation or improper use. Where greater accuracy is required, the following procedures should be observed:

- use modern instruments (manufactured within three years) which are fully checked;
- use certified calibration masses before each weighing operation;
- make at least two measurements;
- consider the use of two instruments installed together and make weighings on both units, otherwise use calibrated masses with a certificate of calibration from a recognised institute.

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WARNING! If the instrument STOPS WORKING because of incorrect operations on the keyboard, DISCONNECT THE MAIN PLUG, then RECONNECT AGAIN pressing simultaneously ON/OFF key [2].

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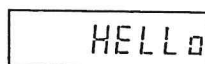
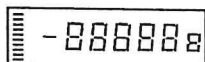
A rectangular digital display showing the word "HELLO" in a simple, blocky font.

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A rectangular digital display showing a test pattern consisting of a series of dashes and numbers, specifically "- 888888 8".

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A rectangular digital display showing the number "0.0" in a simple, blocky font.

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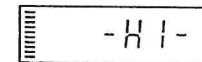
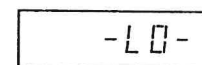
A rectangular digital display showing the message "-LO-" in a simple, blocky font.

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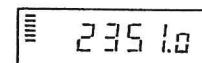
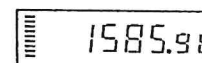
A rectangular digital display showing the number "2351.0" in a simple, blocky font.

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A rectangular digital display showing the number "0.00" in a simple, blocky font.

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A rectangular digital display showing the number "1585.98" in a simple, blocky font.

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This card is of no value if not properly filled-in and it shall be shown to take advantage of the warranty.
It confers either "free-of-charge" replacement or repair of the components recognized faulty from the origin, provided the instrument is returned "freight-prepaid" to the distributor.

The following are excluded:

- damages caused during the transit
- faulty power connection
- accidental unforeseen faults
- damages caused by improper use and/or not in conformity with the enclosed instruction
- glass parts.

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