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***Before using the equipment read
this user guide***

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Even if this device is provided with protections, the use of equipment without following the described procedures could produce a wrong working and equipment's damages.

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This equipment has been designed, built and tested following the European Directive 89/336 CEE (EMC electro-magnetic compatibility), are in conformity with the European Directive 73/23/CEE (BT low tension), and the normative CEI EN 61010-1

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TABLE OF CONTENTS

1	DESCRIPTION	page 11
2	OPERATING FEATURES	page 11
3	INSTALLATION	page 12
4	INSTRUCTION FOR USE	page 12
5	MAINTENANCE	page 12
6	WORKING DEFECTS	page 12
7	GUARANTEE	page 13
8	DOCUMENT VALIDITY	page 13
9	SPARE PARTS	page 13
10	WIRING DIAGRAM	page 13
11	PARTS LIST	page 14
12	DECLARATION OF CONFORMITY	page 15

1 - DESCRIPTION

The Heating Magnetic Stirrer mod. 710/R has a top plate capable to carry high capacity vessels, it is particularly suitable for continuous operations where heating is required. Electronic regulation of magnet speed from 50 to 1000 revolutions per minute. Continuous and noiseless operation. Maximum resistance regulation: 330 W. Outer casing in stainless steel.

2 - OPERATING FEATURES

2.1 - TECHNICAL SPECIFICATIONS

- ◆ Overall dimensions: 195 (width) x 210 (depth) x 125 (height) mm.
- ◆ Working surface dimensions: 160 x 160 mm.
- ◆ Weight: 3 Kg

2.2 - ELECTRICAL SPECIFICATIONS

- ◆ Power supply: 230 V - 50 Hz.
- ◆ Electrical input: 330 W
- ◆ Overload fuses: 2 fuses x 2 A F (5x20) mm.
- ◆ Connection outlet: 10 A

2.3 - CLIMATIC CONDITION FOR STORAGE, PACKING AND USE

- ◆ Altitude up to 2000 m
- ◆ Temperature 5 °C to 40 °C
- ◆ Maximum relative humidity 80% for temperatures up to 31 °C decreasing linearly to 50% relative humidity at 40 °C
- ◆ Mains supply voltage fluctuations not to exceed $\pm 10\%$ of nominal voltage
- ◆ Transient overvoltages II
- ◆ Pollution degree 2

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3 - INSTALLATION

In order to have a functional working position, it is necessary to put the item on a stable support, far from possible obstacle that can influence the normal working area. To move it lift from lower part. Don't use in dangerous atmosphere. Don't use to stir dangerous material or toxic gases that can bring about implosion or explosion, without precaution.

Before the connection of the equipment check that:

- ◆ The rated voltage value indicated on the equipment nameplate corresponds to the actual mains values.
- ◆ The power supply has an effective connection on ground.
- ◆ The electric controls are in a not-working position.
- ◆ Plug in the power supply cable.

4 - INSTRUCTION FOR USE

Turn the main switch on, green led on (Pos 1, Fig.1).

Turn the speed knob till the max. value, put it on the desired speed value turning it in a non-clockwise direction (Pos. 2, Fig. 1).

In order to regulate the temperature use the appropriate knob (Pos.3, Fig.1).

5 - MAINTENANCE

Cleaning: box powdered painted, use ammonia's products.

To remove contaminated substance use sodium hypochlorite (5%)

The maintenance has to be carried out by experts.

To look at the good state of the item and of the plug.

In order to avoid any contact with hot parts, the operations have to be done in total safety conditions (disconnect the plug from the supply power cable).

6 - WORKING DEFECTS

In case of lack of tension, item stop working.

In case of failure to ignite:

- ◆ check the fuses
- ◆ check the outlet is in a good state

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7 - GUARANTEE

- ◆ The good working of the described product is guaranteed from ASAL for a period of 12 months from the purchase date.
- ◆ The guarantee includes the repair or the free exchange of the instrument components that ASAL technicians admit to be defective in terms of manufacture or material (carriage and packing costs to be paid by the buyer).
- ◆ The guarantee will be invalidated by negligence, wrong installation and uses which are not observed in this instructions book, incident and tampering or transfer damage (excluding when it is made by our authorized people).
- ◆ The guarantee extension is no longer valid after repair.
- ◆ In case of restitution, the equipment has to be returned clean in order to avoid any problem for our personnel. We will refuse to draw back all the instruments that do not observe these sanitary regulations.
- ◆ Fuses are not included in the guarantee.

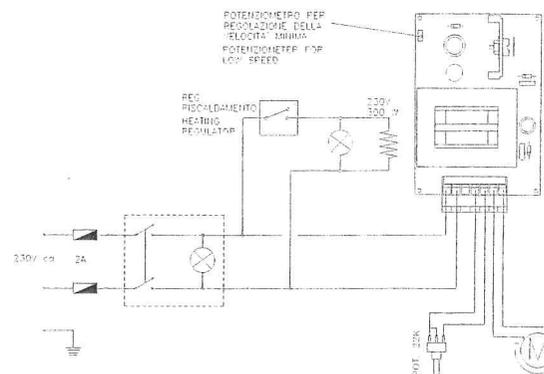
8 - DOCUMENT VALIDITY

ASAL reserves the right to review this document without any notice (any variations do not prejudice the observation of the directives).

9 - SPARE PARTS

For any spare part needed, contact the builder (see address and telephone number on the last page).

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