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Declaration of compliance

We **KARTELL spa**
address **Via Delle Industrie, 1**
20082 NOVIGLIO (MI)
Italy

Under our responsibility declare that the product conforms to inforce prescriptions on safety and hygiene of machines directive 89/392/CEE, low voltage directive 73/23/CEE, electromagnetic compatibility directive 89/336/CEE plus modifications and that the documents listed in the annex V are available at Velp's offices as per the machine directive.

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Instruction for use: vibrating shaker type TK3S

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IMPORTANT:

READ THE INFORMATION CONTAINED IN THE PRESENT MANUAL BEFORE USING THE UNIT.
THE MANUFACTURER DOES NOT ACKNOWLEDGE ANY RESPONSABILITY FOR AN IMPROPER USE OF THE EQUIPMENT, NOT RESPONDING TO DIRECTIONS FOR USE.

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The labels applied to the unit warn the user on the dangers to which he is exposed during the use or the maintenance. The labels must be left on the unit and substituted if they are no more readable.

Warn of danger

To prevent accidents which could be caused by the unit

Start the operation

Do not start to use the unit before reading directions for use and maintenance

SAFETY RULES

- 1) *Hold correctly the test tube and lean it on the rubber cup which rotating transmits vibration to the contained material.*
- 2) *If a fuse is to be replaced, first of all unplug the unit.*

CLEANING

Always unplug the unit before cleaning. The heating plate must be cool. Use a damp cloth and not flammable, not aggressive detergents.

PERSONAL PROTECTION EQUIPMENT

The equipment used for personal protection must be compatible with the dangers due to the working materials and from the glass vessel.

MAINTENANCE

According to the law on products guarantee our instrument must be repaired at our factory, save different agreements with our agents.

GUARANTEE

Starts from the date of delivery note and is referred to the number of register of the single unit.

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
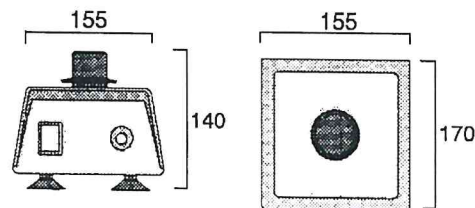
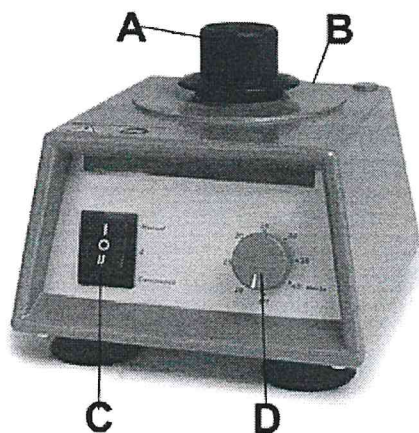
Instruction for use

1. Connect the power cable to mains 230V/50Hz (115V/60Hz) plus ground.


2. When the switch (C) is in "manual" position the unit starts when a slight pressure is exerted on the neoprene small cup (A) and stops by releasing the pressure. When the switch (C) is in "continuous" position the shaking frequency can be selected by turning the electronic regulator knob (D).

3. The apparatus is equipped by 2 fuses connected in the socket (B) located on the rear panel. A blown fuse can be changed after removing rear connection socket and opening the small cover, using the tip of a screwdriver.

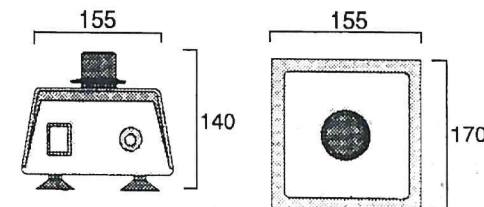4. The equipment is manufactured in accordance with the following safety standards IEC 1010 (1990) + modification 1 (1991).

Power	45 W
Weight	2,15 kg
Supply	230V/50 Hz
Supply	115V/60 Hz

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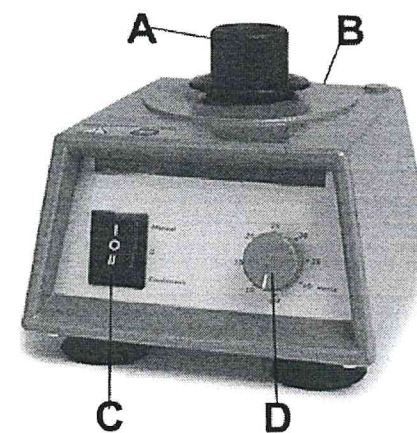Electric scheme 230V/50Hz (115V/60Hz)

- 1 5x20mm retarded fuse 1 A (3.15A)
- 2 Electric motor
- 3 Speed controlling electronic card
- 4 Potentiometer
- 5 Double polar deviator

Note: